

## HybriSlip Hybridization Covers

### INSTRUCTIONS FOR USE

1. With forceps remove the scored polyethylene liner to expose a HybriSlip.
2. Remove the HybriSlip from the unscored liner and apply to specimen as you would a conventional glass cover slip.

---

HybriSlips are RNase free and should be handled with forceps to prevent contamination.

HybriSlips are intended to be used for incubation steps of *in situ* assays and are chemically resistant to the reagents commonly used.

HybriSlips do not require silanization to prevent probes from sticking. Either surface may be used.

